

Contrast RED Solution

for Immunohistochemistry

Catalog No.
71-00-05

Size
50 mL

DESCRIPTION

Contrast RED Solution for Immunohistochemistry contains a nuclear fast red stain. It is designed for use as a counterstain with the HistoMark immunohistochemical staining systems.

FORM/STORAGE/STABILITY

Contrast RED is a 50 mL solution. Store tightly capped, refrigerated at 2 - 8°C. Stable for at least one year stored at 2 - 8°C.

USE

Product is provided at working concentration; no dilution is required.

After immunoassay is complete and tissue sections or other cellular preparations have been reacted with substrate solution, rinse slide with running tap water for 2 - 3 minutes. Counterstain specimen with Contrast RED Solution for 5 - 10 minutes.

After incubation with Contrast RED, rinse specimens